DISCOVERY

Comfort

ENGLISH	1 – 16
DEUTSCH	17 – 34
FRANÇAIS	35 – 52
ESPAÑOL	53 – 68
POLSKI	69 – 84
РУССКИЙ	85 – 100




CE **IVD**

HTL
HTL LAB SOLUTIONS

PZ HTL S.A. is in possession of a registered and certified quality management system which includes the development, production, and sales of high quality Liquid Handling products.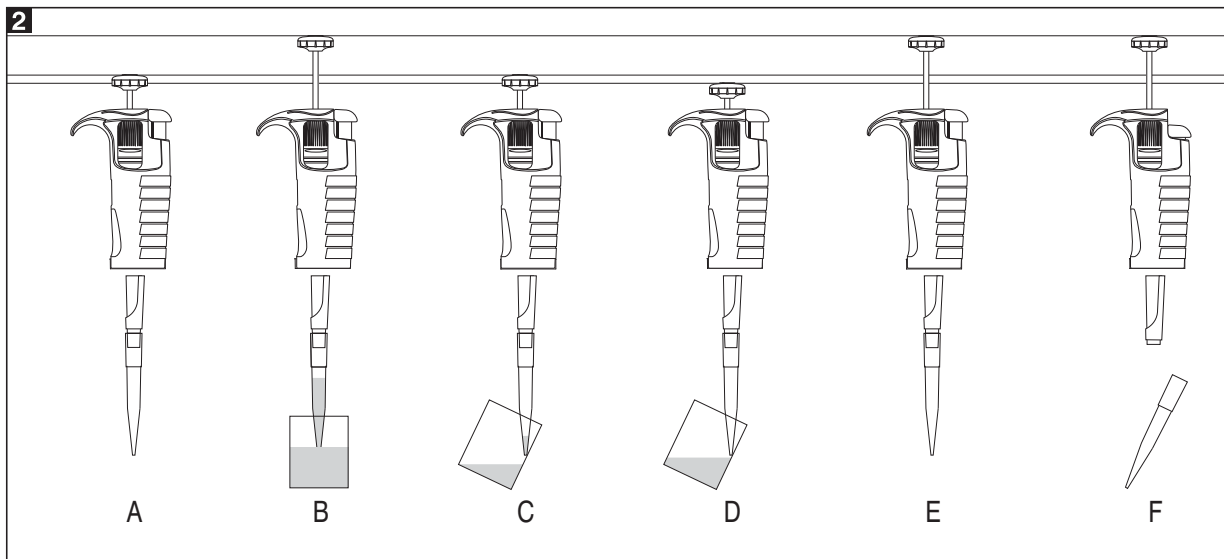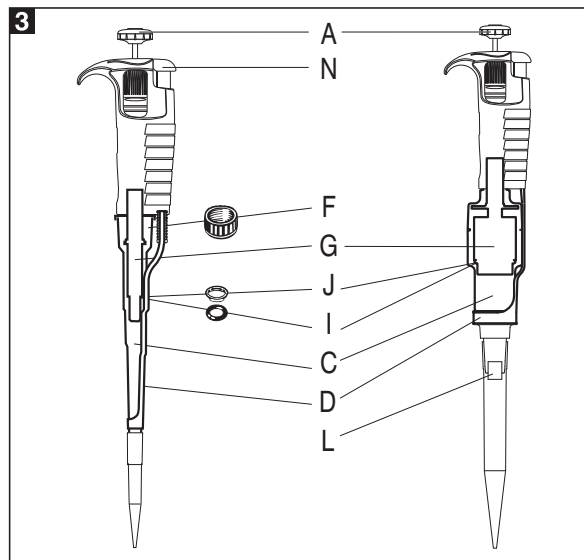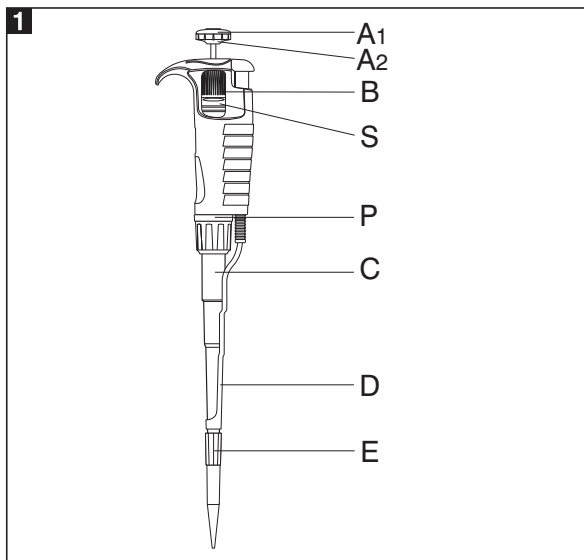

CONTENTS

1 - INTRODUCTION
2 - SETTING THE VOLUME
3 - ASPIRATING AND DISPENSING LIQUIDS
4 - PRE-RINSING
5 - DENSE AND VISCOUS LIQUIDS
6 - HTL PIPETTE TIPS
7 - RECOMMENDATIONS
8 - RECALIBRATION
9 - TROUBLESHOOTING
10 - CLEANING AND STERILIZATION
11 - PIPETTE KIT
12 - SPARE PARTS

1 - INTRODUCTION

The **DISCOVERY Comfort** is a volumetric instrument designed to measure and transfer liquids precisely and safely. It can measure and transfer volumes from 0.1 μl to 10000 μl depending on the model.

The **DISCOVERY Comfort** pipettes are available in two versions which differ in the shaft color:

DV - the shafts are made in a light grey colour for all volumes of pipettes.

D - non standard version on special request only. The shafts are made in red, yellow, green, blue, and white colour depending on the volume range of the pipette.

The **DISCOVERY Comfort** pipettes are equipped with a digital counter which shows the pipetting volume. The set volume is visible in the window on the handle. The setting of the volume is done by turning of the pipetting pushbutton knob (Fig. 1A2) or by turning the adjustment knob (Fig.1B). The pipette design allows the user to lock the volume setting by pushing the locking ring upwards (Fig.1S). The position of the ring is indicated by the symbols located on the handle. The range of the volume of the aspired liquid is shown on the pipetting pushbutton (Fig. 1A1).

The **DISCOVERY Comfort** pipettes are produced in 10 models, covering the volume range from 0.1 μl to 10000 μl .

Model	Volume range [μl]
DV2, D2	0.1 - 2
DV10, D10	0.5 - 10
DV20, D20	2 - 20
DV50, D50	5 - 50
DV100, D100	10 - 100
DV200, D200	20 - 200
DV250, D250	50 - 250
DV1000, D1000	100 - 1000
DV5000, D5000	1000 - 5000
DV10000, D10000	1000 - 10000

DV2, DV10, D2, D10	Measurement and transfer of micro-volumes, DNA sequencing and enzyme-assay applications.
DV20, DV50, DV100, DV200, DV250, DV1000, D20, D50, D100, D200, D250, D1000	Measurement and transfer of general aqueous solution, acids and bases
DV5000, DV10000, D5000, D10000	Measurement and transfer of large volumes.

The liquid is aspirated into disposable tips attached to the pipette shaft. Disposable tips ensure maximum safety and eliminate possibility of crosscontamination of the liquid samples.

To protect the user from contaminated tips the **DISCOVERY Comfort** is equipped with a built-in tip ejector, (Fig. 1D).

The construction of the ejector enables the user to set up the length. The adjustable tip ejector accommodates every type of tips available on the market. When using narrow tubes, it may be necessary to remove the tip ejector. It is simply removed by pulling down.

Adjusting of the Tip Ejector Length

- in 2-1000 μl pipettes (Fig. 6A).

The regulated "H" spacers, included in the box allow for regulating the length of tip ejector by +1mm or +2mm. An "H0" spacer is inserted on a standard basis. The outside shape of the spacer identifies the adjustment.

- in 5000 and 10000 μl pipettes (Fig. 6B).

The length of tip ejector is regulated by screwing in or screwing out the tip ejector stem with a screwdriver. Turn the screwdriver counter clockwise to increase the length of tip ejector, and clockwise to reduce the length of tip ejector. The ejector can be adjusted in the range of 5 mm.

If above described method of ejector adjustment is not sufficient or the diameter of the ejector opening is too large to eject the tip it is necessary to put the ejector cap "M" onto the ejector, (Fig. 6C).

- in 2 and 10 μl pipettes (Fig. 6D)

The 2 μl - 10 μl pipettes may require the use of an ejector cap "M" to efficiently eject certain brands of pipette tips. Simply place the cap, supplied with the pipette, on the bottom of the pipette shaft and slide the cap upwards until it surrounds the bottom of the tip ejector.

The **DISCOVERY Comfort** is a high quality instrument which offers excellent accuracy and precision.

The accuracy and precision (repeatability) of liquid volume depend on the quality of tips used. The values for accuracy and precision given in the table below were obtained using **HTL** tips. Only those tips guarantee correct operation of the pipettes and ensure accuracy and precision of liquid sampling.

Model	Cat. no.	Volume [μl]	Accuracy [%]	Precision [%]	Fit to tips μl	Shafts color for D
DV2 D2	4041 4051	0.2 1.0 Max 2.0	± 12.0 ± 2.7 ± 1.5	± 6.0 ± 1.3 ± 0.7	10	red
DV10 D10	4042 4052	Min 0.5 5.0 Max 10.0	± 4.0 ± 1.0 ± 0.5	± 2.8 ± 0.6 ± 0.4		
DV20 D20	4043 4053	Min 2 10 Max 20	± 3.0 ± 1.0 ± 0.8	± 1.5 ± 0.5 ± 0.3	200	yellow
DV50 D50	4047 4057	Min 5 25 Max 50	± 2.5 ± 1.0 ± 0.8	± 2.0 ± 0.6 ± 0.4		
DV100 D100	4044 4054	Min 10 50 Max 100	± 1.6 ± 0.8 ± 0.8	± 0.80 ± 0.24 ± 0.20		
DV200 D200	4045 4055	Min 20 100 Max 200	± 1.2 ± 0.8 ± 0.6	± 0.60 ± 0.25 ± 0.20		
DV250 D250	4040 4050	Min 50 125 Max 250	± 1.0 ± 0.8 ± 0.6	± 0.4 ± 0.3 ± 0.3	300	green
DV1000 D1000	4046 4056	Min 100 500 Max 1000	± 1.6 ± 0.7 ± 0.6	± 0.40 ± 0.20 ± 0.15	1000	blue
DV5000 D5000	4048 4058	Min 1000 2500 Max 5000	± 0.6 ± 0.6 ± 0.5	± 0.25 ± 0.20 ± 0.15	5000	white
DV10000 D10000	4049 4059	Min 1000 5000 Max 10000	± 2.5 ± 0.8 ± 0.5	± 0.6 ± 0.3 ± 0.2	10000	

The accuracy and precision are obtained with **HTL** tips, using a gravimetric method, performing at least 10 measurements of distilled water at a temperature of $20 \pm 1^\circ\text{C}$, according to EN ISO 8655 standard.

The pipette design enables the user to perform the recalibration process according to the rules presented in section 8.

2 - SETTING THE VOLUME

The volume display shown by the counter has three digits, which should be read from top to bottom. In addition the lowest counter drum is printed in a scale which enables the setting of the volume in the minimum graduation range. Examples of the meanings of the black and red digits:

Pipettes **DV2, D2**

Red figures at the bottom = $1/100 \mu\text{l}$
Increment = $0.002 \mu\text{l}$

	DV2, D2
	1
red	2
red	5
	1.25 μl

Pipettes **DV10, DV20, D10, D20**

Red figures at the bottom = $1/10 \mu\text{l}$
Increment = $0.02 \mu\text{l}$

	DV10, D10	DV20, D20
	0	1
	7	2
	5	5
	7.5 μl	12.50 μl

Pipettes **DV50, DV100, DV200, DV250, D50, D100, D200, D250**

Black figures only = $1 \mu\text{l}$
Increment = $0.2 \mu\text{l}$

	DV50, DV100 D50, D100	DV200, D200 DV250, D250
	0	1
	7	2
	5	5
	75 μl	125 μl

Pipettes **DV1000, DV5000, D1000, D5000**

Red figures at the top - ml
Increment =

	DV1000 D1000	DV5000 D5000
	0	1
red	7	2
red	5	5
	0.75 ml	1.25 ml

2 μl 10 μl

Pipettes **DV10000, D10000**

Red figures at the top - ml
Increment = $20 \mu\text{l}$

	DV10000, D10000
	0
red	7
red	5
	7.5 ml

The volume of the pipette is set by the knob in the pipetting pushbutton (Fig. 1A2) or by the black adjustment knob (Fig. 1B). The volume adjustment can be performed when the locking ring is set in the lower position (Fig.5A).

When the desired volume is selected, the locking ring should be set in the upper position (Fig.5B). To attain the maximum accuracy, set volume must be approached from a higher value by diminishing counter readings.

- If the desired volume is lower than set volume shown by the counter, the operator should turn the pipetting pushbutton (Fig. 1A2) or the black adjustment knob (Fig. 1B) to the direction diminishing counter readings to the required volume. Before achieving the required volume slowly rotate the knob and observe carefully diminishing reading to avoid accidentally passing the setting value.
- If the desired volume is higher than set volume shown by the counter, the operator should turn the pipetting pushbutton (Fig. 1A2) or the black adjustment knob (Fig. 1B) increasing the value until the lower figure wheel comes 1/3 of a turn beyond the required setting and then slowly backward until the setting reaches the desired volume. Make sure not to pass the setting value.

If the knob is accidentally turned too far, the process must be repeated. The desired volume must always be set from the higher value in the order of decreasing value.

Following volume adjustment, set the locking ring into the upper position, thus locking the knob and preventing accidental volume change.

3 - ASPIRATING AND DISPENSING LIQUIDS

Place a tip on the shaft of the pipette. See Section 6 for the appropriate tip. Press the tip on firmly using a slight twisting motion. This will ensure an airtight seal.

Important: Never aspirate liquids into the DISCOVERY Comfort without a tip attached.

Aspiration

Press the pushbutton to the first positive stop, (Fig. 2A). Holding the pipette vertically, immerse the tip into the sample liquid. The depth to which the tip is immersed in the sample liquid depends on the model.

Model	Immersion depth (mm)
DV2, D2	≤ 1
DV10, D10	≤ 1
DV20, DV50, DV100, D20, D50, D100	2 ÷ 3
DV200, DV250, DV1000, D200, D250, D1000	2 ÷ 4
DV5000, D5000	3 ÷ 6
DV10000, D10000	5 ÷ 7

Release the pushbutton slowly and smoothly to aspirate the sample, (Fig. 2B). Wait one second and then withdraw the tip from the liquid. When the pipette tip is immersed not as deeply as the recommended depth or when the pipetting pushbutton is rapidly released air may enter the disposable tip.

Avoid touching the orifice of the tip.

Dispensing

- Place the end of the tip against the inside wall of the vessel at an angle of 10 to 40 degrees.
- Press the pushbutton smoothly to the first stop, (Fig. 2C). Wait one second.
- Press the pushbutton to the second stop to expel any remaining liquid, (Fig. 2D).
- Keeping the pushbutton depressed to the very end, remove the pipette by drawing the tip against the inside surface of the receiving vessel.
- Release the pushbutton to its starting position, (Fig. 2E).
- Eject the tip by pressing the tip ejector button, (Fig. 2F). Remember to change the tip whenever a different kind of liquid is to be sampled.

Filters

A replaceable filter installed in a seat in the bottom part of the shaft is used in 5000 µl and 10000 µl pipettes (Fig. 3L). The filter prevents the aspirated liquid from entering into the shaft and thus from polluting the inside of the shaft and the piston. Using the filter is especially important when aspirating and dispensing large volumes of liquid.

If the filter becomes wet during liquid aspiration it should be replaced with a new one.

4 - PRE-RINSING

When pipetting liquids of higher viscosity or lower surface tension than water (e.g. sera or organic solvents),

a film of liquid is formed on the inside wall of the pipette tip. This film can create an error. Since the film remains relatively constant in successive pipetting operations with the same tip, this error can be avoided by forming the film before transferring the first sample. This is done by aspirating a sample and dispensing it back into the same vessel. Since the film is already formed, all of the following samples will have better accuracy and repeatability.

This pre-rinsing operation should be repeated when the volume to be aspirated is changed or when a new tip is used.

5 - DENSE AND VISCOUS LIQUIDS

The **DISCOVERY Comfort** specifications of accuracy and precision are based on pipetting distilled water. The handling of liquids with physical qualities of density, viscosity and surface tension differing extremely from water may need a gravimetrically checked compensation of the volume setting. Normally the degree of error resulting from heavy or viscous liquids is negligible if the pipetting is done slowly and carefully. It is most important to give the liquids some time to follow the change of pressure by holding the pipette tip in its position for at least 2 sec. after the aspiration and the blow out stroke.

If in extreme cases this method of operation does not result in accurate values, a compensation could be achieved as follows:

Weigh the liquid pipetted when the **DISCOVERY Comfort** is set to the nominal value. Then calculate the set-off from the nominal value:

$$\text{Corr: val.} = 2 \times \text{nom. val.} - \frac{m}{\gamma}$$

m - weight of the sample

γ - density of liquid

Check this operation once again and correct if necessary. Note the corrected value for further pipetting the same kind of liquid.

6 - HTL PIPETTE TIPS

HTL tips are made from high performance polypropylene and their quality guarantees the precision and accuracy associated with the **DISCOVERY Comfort**. Strict control is maintained throughout the manufacturing process to ensure the highest quality.

The accuracy and precision figures for the **DISCOVERY Comfort** are only guaranteed when **HTL** tips are used. The use of inferior quality tips will seriously degrade the performance of the **DISCOVERY Comfort**.

Tips 10

These tips are used for volumes between 0.1 μl and 10 μl . They are used with the DV2, DV10, D2 and the D10 models, which are equipped with the red pushbutton.

Tips 200

These tips are used for volumes between 2 μl and 200 μl . They are used with the DV20, DV50, DV100, DV200, D20, D50, D100 and the D200 models, which are equipped with the yellow pushbutton.

Tips 300

These tips are used for volumes between 50 μl and 250 μl . They are used with the DV250 and the D250 models, which are equipped with the green pushbutton.

Tips 1000

These tips are used for volumes between 100 μl and 1000 μl . They are used with the DV1000 and the D1000 models, which are equipped with the blue pushbutton.

Tips 5000

These tips are used for volumes between 1000 μl and 5000 μl . They are used with the DV5000 and the D5000 models, which are equipped with the white pushbutton.

Tips 10000

These tips are used for volumes between 1000 and 10000 μl . They are used with the DV10000 and the D10000 models, which are equipped with the white pushbutton.

7 - RECOMMENDATIONS

Observing the following recommendations will ensure maximum possible accuracy and precision of liquid sampling.

- When pipetting the **DISCOVERY Comfort** should have the volume setting locked with the locking ring in the upper position.
- Make sure to operate the **DISCOVERY Comfort** slowly and smoothly.
- The depth of immersion in the sample liquid should be the minimum necessary and should remain constant during aspiration.

- The **DISCOVERY Comfort** should be held in a vertical position.
- Change the tip when volume setting is changed or when a different liquid is to be aspirated.
- Change the tip if a droplet remains on the end of the tip from the previous pipetting operation.
- Each new tip should be pre-rinsed with the liquid to be pipetted.
- Liquid should never enter the **DISCOVERY Comfort** shaft. To prevent this:
 - Press and release the pushbutton slowly and smoothly.
 - Never turn the pipette upside down.
 - Never lay the pipette on its side when there is liquid in the tip.
- Never force the volume setting beyond its recommended limits.
- When pipetting liquids with temperatures different from the ambient temperature, it is recommended to pre-rinse the tip several times before use.
- Do not pipette liquids with temperatures above 70°C.
- When pipetting acids or corrosive solutions which emit vapours, it is recommended to disassemble the shaft and to rinse the piston and seal with distilled water after finishing the pipetting operation.

8 - RECALIBRATION

HTL pipettes are calibrated by gravimetric method, using **HTL** tips and distilled water, at the temperature $20 \pm 1^\circ\text{C}$, according to EN ISO 8655 standard.

If during pipette operation you find that the accuracy error (the difference between the real aspirated volume and the preset volume) exceeds the permissible value given in the table in section 1, the pipette recalibration procedure should be carried out.

Before starting the recalibration it is necessary to check whether the following requirements have been fulfilled during error determination:

- the ambient temperature, and the temperature of the pipette, tips and water was identical
- the density of the liquid used is close to that of distilled water
- a balance with appropriate sensitivity has been used.

Volume checked [μl]	Balance sensitivity [mg]
0.1 - 10	≤ 0.001
10 - 100	≤ 0.01
> 100	≤ 0.1

- mg/ μl conversion factor has been taken into account
- the requirements given in sections 3 and 7 have been fulfilled

If the above conditions are satisfied and the accuracy error for selected volume given in section 1 exceeds the permissible value, the pipette recalibration procedure should be carried out.

The recalibration can be performed within one full turn of the key to the right or to the left only.

Recalibration conditions:

- Ambient temperature and the temperature of the pipette, tips and liquid should be within the range $20-25^\circ\text{C}$ and stabilised during weighing within $\pm 0.5^\circ\text{C}$
- Measurements should be conducted using distilled water
- Balance sensitivity should be suitable for the volume to be controlled

Recalibration procedure:

- Set the dose volume depending on the pipette volume according to the following table:

Model	Range of the pipette volumes [μl]	Preset volume [μl]	Permissible volumes [μl]	Volume change ΔV for full turn of the calibration key [μl] (24 increments)
DV2, D2	0.1 - 2	0.2	0.176 - 0.224	0.06
DV10, D10	0.5 - 10	0.5	0.48 - 0.52	0.33
DV20, D20	2 - 20	2	1.94 - 2.06	0.63
DV50, D50	5 - 50	5	4.875 - 5.125	2.50
DV100, D100	10 - 100	10	9.84 - 10.16	2.50
DV200, D200	20 - 200	20	19.76 - 20.24	6.30
DV250, D250	50 - 250	50	49.5 - 50.5	6.30
DV1000, D1000	100 - 1000	100	98.4 - 101.6	25.00
DV5000, D5000	1000 - 5000	1000	994 - 1006	125.00
DV10000, D10000	1000 - 10000	1000	975 - 1025	250.00

- Perform 5 aspirations, weigh each one and calculate the average value of the aspirations

- Calculate average aspirated volume in μl multiplying the average aspiration amount [mg] by the distilled water density coefficient [$\mu\text{l}/\text{mg}$], which depends on temperature and pressure according to the following table:

Temperature [°C]	Pressure [kPa]		
	95.0	101.3	105.0
20	1.0028	1.0029	1.0029
21	1.0030	1.0031	1.0031
22	1.0032	1.0033	1.0033
23	1.0034	1.0035	1.0036
24	1.0037	1.0038	1.0038
25	1.0039	1.0040	1.0040

If the average aspirated volume exceeds the permissible value, the following should be done:

- Remove the pipetting pushbutton, (Fig. 4A),
Warning: The pipetting pushbutton consists of 2 parts: the knob (Fig. 1A2) and the pushbutton (Fig. 1A1). After removal of the pushbutton, both parts are separated.
- Holding the volume setting knob to protect it against rotation, insert the calibration key into the cuts of the calibration screw, (Fig. 4B),
- Turn the key clockwise to reduce the aspirated volume, or counter-clockwise to increase the volume, (Fig. 4C). One full turn of the calibration key changes the pipette aspiration volume by the amount given in the table.
- Take out the key and fix the pipetting pushbutton (Fig. 4D). The pipetting pushbutton should be fixed by placing the knob on the arbor first (Fig. 1A2) and then the pushbutton (Fig. 1A1).

Determine the average aspirated volume. The average volume should be within the permissible range given in the table. If the volume exceeds the values stated, the recalibration procedure should be repeated.

When pipetting liquids with physical properties considerably different from those of water, follow the rules given in section 5.

More information on calibration procedure can be found on www.htl.com.pl

9 - TROUBLESHOOTING

If you notice an improper pipette operation identify the cause and eliminate the fault. To do this, follow the instruction in the sequence provided. Replacement of parts should be required only occasionally, and should not occur under normal pipette use.

Droplets of liquid remain in the pipette tip.

- The tip is emptied too fast.
Decrease the speed of pressing the pipette pushbutton.
- The tip wettability has increased due to extensive use.
Replace the tip with a new one.

Droplets of air appear in the liquid aspirated into the tip.

- The pipette tip immersion is too shallow.
Immerse the tip deeper according to the instructions.
- The pipette tip is incorrectly pressed onto the pipette shaft.
Press the pipette firmly.
- The tip is damaged or worn out due to extensive use.
Replace the tip with a new one.

The pipette incorrectly aspirates the liquid or liquid drops out from the tip.

- The pipette tip is incorrectly pressed onto the pipette shaft.
Press the pipette tip firmly.
- The shaft nut is loose (Fig. 3F) in the models DV/D2-DV/D1000
Tighten the shaft nut.
- The sealing surface of the shaft is cracked or scored.
Remove the tip ejector. Unscrew the shaft nut, inspect the shaft and the piston assembly. Replace the damaged parts (see Section 12). When reassembling the pipette, the nut should be hand tightened. In the models DV/D2, DV/D10 and DV/D20, the damage of the shaft may also cause a damage of the piston assembly. Replace the damaged parts (see Section 12). When reassembling the pipette, the nut should be hand tightened.
To remove the tip ejector in models DV/D5000 and DV/D10000, remove the ejector pushbutton (Fig. 3N) and using a screwdriver unscrew the tip ejector by turning the screwdriver counter-clockwise.

- Damage to the piston or seal due to prolonged use with the aggressive liquids.

Disassemble the pipette as described above. Replace the piston, seal and O-ring (see Section 12). Rinse the inside of the shaft in distilled water and dry. Lubricate the seal and O-ring with the lubricant, that has been included with each pipette.

The replacement of the piston requires conducting of calibration procedure.

Note: The parts of DV2/D2 and DV/D10 pipette should be lubricated evenly with a minimum amount of lubricant.

- The pipette is reassembled improperly.
Disassemble the pipette and reassemble it, observing the proper sequence of steps (Fig. 3).
- No lubricant on the sealing elements.

Remove the tip ejector. Unscrew the shaft nut, remove the shaft, piston assembly, seal and O-ring. Rinse the removed parts in distilled water and dry thoroughly. Lightly lubricate the inside surfaces of the seal and the O-ring with the included lubricant. Reassemble the pipette in the reverse order.

- Contamination of the inside of the pipette caused by extensive aspiration of chemically aggressive liquids or because liquid got inside the pipette.

Remove the tip ejector. Unscrew the nut, remove the shaft, piston assembly, seal and O-ring. Rinse the removed parts with distilled water and dry thoroughly. Lightly lubricate the inside surfaces of the seal and the O-ring with the lubricant. Reassemble the pipette in the reverse order.

If you find an increase in the pipetting force, which could happen after repetitive autoclaving of the pipette: Remove the tip ejector. Unscrew the shaft nut, and then remove the shaft, piston assembly, seal and O-ring. Rinse the removed parts in distilled water and dry. Lubricate the internal surfaces of the seal and O-ring with lubricant that has been included with each pipette. Reassemble the pipette in opposite order.

Note: All parts of the pipette can be autoclaved (see Section 10)

The shaft of the 5000 and 10000 models should be autoclaved without the filter.

If the problem continues after carrying out the above steps, contact your **HTL** representative.

Before returning the pipette, please ensure that the pipette is completely free of any chemical, radioactive or microbiological contamination which could pose a threat during transport and repair. As far as it is possible, clean the pipette.

10 - CLEANING AND STERILIZATION

Cleaning

External surfaces of the pipetting pushbutton, the ejector pushbutton, the handgrip, the shaft nut and the adjustment knob may be cleaned using a cloth dampened in isopropyl alcohol. The remaining parts removed from the pipette during pipette disassembly may be washed with distilled water or isopropyl alcohol.

Sterilization:

The pipette can be sterilized in the autoclave at 121°C for 20 minutes. After sterilization, the pipette should be dried and cooled to room temperature.

It is recommended:

- to sterilize the pipettes in autoclave with an initial vacuum and drying cycle,
- prior to sterilization unscrew the shaft nut slightly in the DV/D2-DV/D1000 pipettes, and unscrew the shaft slightly in the DV/D5000 and DV/D10000. After autoclaving these parts should be screwed tight again.
- to set the locking ring in lower (unlocked) position prior to sterilization

The precision of the results should not alter if the pipetting process and autoclaving are carried out as described in this manual. Because a slight change in the accuracy of the dosage may occur, it is recommended to:

- check the calibration of the pipette after the initial first, third and fifth autoclaving cycles and then after every 10 autoclaving cycles.

11 - PIPETTE KIT

The pipettes are delivered in the kits including:

- Pipette
- Instruction manual
- Calibration key
- Pipette stand
- Ejector regulation spacers
(for pipette models DV/D2 - DV/D1000)
- Ejector cap (for pipette models DV/D2, DV/D10, DV/D5000, DV/D10000)
- Tips
- Identification labels
- Filters (for pipette models DV/D5000, DV/D10000)
- Lubricant
- Color Identification rings

The stand assembly diagram is shown in Fig. 7.

12 - SPARE PARTS

The spare parts indicated in Fig. 3, 4, 5 and 6 that is:

- A: Pipetting pushbutton A1: Pushbutton A2: Knob**
B: Adjustment knob
C: Shaft
D: Ejector
F: Shaft nut
G: Piston assembly
H: Spacer
I: O-ring
J: Seal
K: Calibration key
L: Filter
M: Ejector cap
N: Ejector pushbutton
P: Identification ring
S: Locking ring

can be ordered from your **HTL** representative (type of pipette and name of the part for this pipette should be specified).

Warning: The replacement of the piston requires conducting of calibration procedure according to section 8.

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